

Costar Minicentrifuge. QuickGuide

Warnings The following precautions should be taken when operating or working near the minicentrifuge:

NEVER use the centrifuge in any manner not specified in these instructions.

NEVER operate the centrifuge without a rotor properly attached to the shaft.

NEVER fill tubes while they are in the rotor. Liquid spillage may harm unit.

NEVER put hands in the rotor area unless the rotor is completely stopped.

NEVER move the centrifuge while the rotor is spinning. **NEVER** use solvents or flammables near this or other electrical equipment.

NEVER centrifuge flammable, explosive or corrosive materials.

NEVER centrifuge hazardous materials outside of a hood or proper containment facility.

ALWAYS load the rotor symmetrically. Each tube should be counterbalanced by another tube. Weight of individual tubes may not exceed 3.0 g.

ALWAYS locate the centrifuge within easy access to an electrical outlet.

ALWAYS use only microcentrifuge tubes made from plastic and designed to withstand centrifugal forces of at least 2,000 xg.

Operation: